



illumina Proactive Service connectivity guide

Easily connect your NextSeq™ 500 or NextSeq 550 Sequencing System

Minimize unplanned downtime, run failures, and sample loss, and resolve issues faster if and when they arise

Follow these steps to get connected

- From the Home Screen, select **Manage Instrument**
- From the **Manage Instrument** screen, select **System Configuration** followed by **BaseSpace™ Configuration**
- Move past the Analysis tab to the **BaseSpace** tab and check the box next to **Send instrument performance data to Illumina**
- Click **Save**

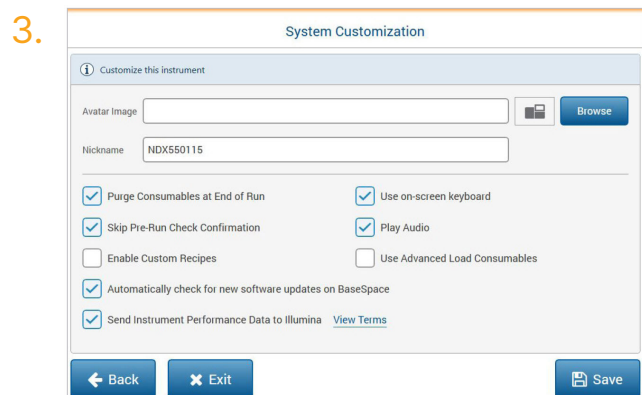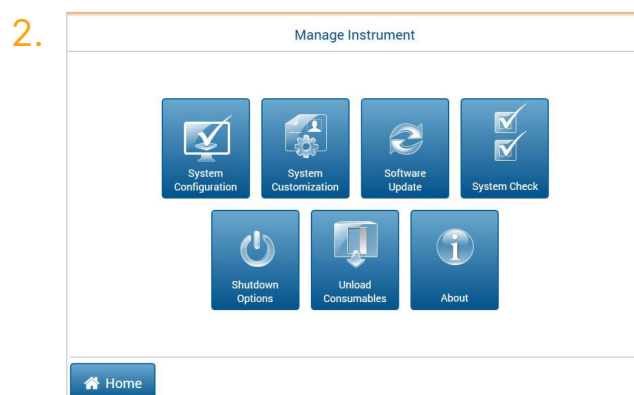

For more information, contact your local field applications scientist or field service engineer

illumina • 1.800.809.4566 toll-free (US) • +1.858.202.4566 tel • techsupport@illumina.com • www.illumina.com

© 2021 Illumina, Inc. All rights reserved. All trademarks are the property of Illumina, Inc. or their respective owners. For specific trademark information, see www.illumina.com/company/legal.html. M-GL-00018